

Brookhaven National Laboratory/ LIGHT SOURCES DIRECTORATE			
<b>Subject:</b>	<b>VACUUM PROCEDURES FOR BEAMLINE X-7A</b>		
<b>Number:</b>	LS-OPS-0133	<b>Revision:</b>	B
		<b>Effective:</b>	08/05/2010
			<b>Page 1 of 3</b>
Prepared By: R. Greene	Reviewed By: J. Klug	Approved By: S. Ehrlich	Approved By: E. Hu

\*Approval signatures on file with master copy.

**The only official copy of this file is the one on-line in the NSLS website. Before using a printed copy, verify that it is the most current version by checking the document effective date on the NSLS website.**

The following procedures must be followed when bleeding up different beamline sections and when returning these sections to operation (refer to Beam Line Layout Drawing):

## **I. FRONT-END (PROCEDURE TO BE PERFORMED BY NSLS VACUUM GROUP ONLY)**

### **A. Bleed-Up**

1. Notify the Coordinator (Beeper 5824).
2. Refer to Front-End Vacuum Procedures (SLS-07.19-13-1).

### **B. Return to Operation**

1. Notify the Coordinator (Beeper 5824).
2. Refer to Front-End Vacuum Procedures (SLS-07.19-13-1).

## **II. SECTION BETWEEN VALVE 1A AND Be WINDOW 1A**

### **A. Bleed-Up**

1. Notify the Coordinator (Beeper 5824).
2. Close and seal Valve 1A and the Front-End High Vacuum Valve.
3. Hook up turbo pump to this section.
4. Coordinator places Yellow Tag on Valve 1A and the Front-End High Vacuum Valve.
5. Slowly bleed-up with boil-off N<sub>2</sub> while Coordinator monitors Front-End pressure.

### **B. Return to Operation**

1. Pump to  $< 2 \times 10^{-9}$  Torr.
2. Notify the Coordinator (Beeper 5824).
3. Prepare for RGA scan.\*
4. Open Valve 1A into Front-End provided pressure  $< 2 \times 10^{-9}$  Torr downstream of valve.
5. Perform RGA scan.\*
6. If RGA scan or pressure reading (if no RGA scan required) is satisfactory, Coordinator removes Yellow Tag from Valve 1A.
7. Remove any unprotected turbo pump from this section or valve off the turbo pump and place a Yellow Tag on the valve.\*\*

## **III. SECTION BETWEEN Be WINDOW 1A AND A1 WINDOW 2A**

### **A. Bleed-Up**

1. Notify the Coordinator (Beeper 5824).
2. Close and seal Valve 1A.
3. Coordinator places Yellow Tag on Valve 1A.
4. Slowly bleed-up with boil off N<sub>2</sub> while Coordinator monitors pressure between Valve 1A and Be Window 1A.

### **B. Return to Operation**

1. Pump to  $< 10^{-3}$  Torr.
2. Notify the Coordinator (Beeper 5824).
3. Coordinator removes Yellow Tag from Valve 1A provided pressure  $< 10^{-3}$  Torr downstream of Be Window 1A.

<b>Subject:</b>	<b>VACUUM PROCEDURES FOR BEAMLINE X-7A</b>					
<b>Number:</b>	LS-OPS-0133	<b>Revision:</b>	B	<b>Effective:</b>	08/05/2010	<b>Page 2 of 3</b>

**\* NSLS POLICY FOR RGA SCANS (24 HOUR NOTICE REQUIRED)**

An RGA scan is required before returning to operation if there is a major change of hardware in the vacuum system, i.e. changing of samples, mirrors, windows, monochromator crystals or gratings, manipulators, detectors, etc., **with the following two exceptions:**

1. After UHV sample chambers have been bled up for replacing components, an RGA scan will not be required if the chamber pressure is returned to  $< 2 \times 10^{-9}$  Torr and the Front End pressure remains  $< 2 \times 10^{-9}$  Torr when vacuum sections upstream of the chamber are opened into the Front End.
2. If any vacuum section upstream of the bled-up section remains at a pressure of  $< 9 \times 10^{-10}$  Torr as read using a hot-filament ion gauge, when the entire beamline is opened into the Front End, and the Front End pressure does not increase, no RGA is required.

**\*\* NSLS TURBO PUMP POLICY**

An unprotected turbo pump is one not separated from the Front End by a beamline valve which automatically closes in the event of a power loss or a pressure increase at the turbo pump. **No unprotected turbo pump can share a contiguous vacuum with the Front End.**

<b>Subject:</b>	<b>VACUUM PROCEDURES FOR BEAMLINE X-7A</b>		
<b>Number:</b>	LS-OPS-0133	<b>Revision:</b>	B
		<b>Effective:</b>	08/05/2010
			<b>Page 3 of 3</b>

<b>Document Review Frequency</b>
<b>3</b> Years

Review signatures on file  
with master copy of  
controlled document

<b>LIGHT SOURCES DIRECTORATE REVISION LOG</b>		
<b>Document Number:</b>	LS-OPS-0133	
<b>Subject:</b>	VACUUM PROCEDURES FOR BEAMLINE X-7A	
<b>Rev</b>	<b>Description</b>	<b>Date</b>
B	Initial release in to the Controlled Document System.	08/05/2010